# QUICK REFERENCE Conjugated EGF

Pub. no. MAN0009580



Revision 1.0

# Detailed protocol is available online at www.lifetechnologies.com/manuals.

WARNING! For every chemical, read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from www.lifetechnologies.com/support.

Label	Cat. no.	Ex/Em maxima	Quantity
Fluorescein	E3478	494/518	20 µg
Alexa Fluor® 488	E13345	495/519	100 µg
pHrodo™ Green	P35375	509/533	20 µg
Alexa Fluor® 555	E35350	555/565	100 µg
Tetramethylrhodamine	E3481	555/580	20 µg
pHrodo <sup>™</sup> Red	P35374	560/585	20 µg
Texas Red®	E3480	595/615	100 µg
Alexa Fluor® 647	E35351	650/665	100 µg
Biotin	E3477	<300/none	20 µg
Unlabeled (mouse)	E3476	NA	100 µg
Unlabeled (recombinant human)	PHG0311	NA	100 µg

For Research Use Only. Not for use in diagnostic procedures.



# Quick Protocol

- 1. Plate cells and keep on ice for 10 minutes.
- Wash cells with cold Live Cell Imaging Solution (Cat. no. A14291DJ) containing 20 mM Glucose and 1% BSA.
- Add EGF conjugate at 2 μg/mL in Live Cell Imaging Solution containing 20 mM Glucose and 1% BSA.
- 4. Incubate at 37°C for 5–20 minutes.
- Wash cells with Live Cell Imaging Solution (optional for pHrodo<sup>™</sup> Red and pHrodo<sup>™</sup> Green EGF conjugates).
- 6. Analyze cells using appropriate excitation/emission maxima given on page 1.

#### Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

BIGLANGE. LET TECHNOLOGIES CORPORTION AND/ORITIS AFFLUATES) DISCLAM ALL WARRANTES WITH RESPECT TO THE DOLLMENT, DPRESSED OR MAREED, MICLIDNEN DUI NOLIMIETO TO THUSE OF MERCIANAULUUT, PITNESS OR A BARTICULAR PURCE, OR NON-AFFRICANAULUUT, TO THIS DOLLMENT, DPRESSED OR MAREED, MICLIDNES DUI NOLIMIETO TO THUSE OF MERCIANAULUUT, PITNESS OR A BARTICULAR PURCE, OR NON-AFFRICANAULUUT, TO THIS DOLLMENT, DPRESSED OR MAREED, MICLIDNES DUI NOLIMIETO TO THUSE OF MERCIANAULUUT, PITNESS OR A BARTICULAR PURCE, OR NON-AFFRICANAULUUT, TO THIS DOLLMENT, DPRESSED OR MAREED, MICLIDNES DUI NOLITOTO TO THUSE OF MERCIANAULUUT, PITNESS OR A BARTICULAR PURCE, OR NON-AFFRICANAULUUT, TOTAL MARANTI, ON UNDERNISTI ON AMARTINI, ON UNDERLES AND AFFRICANAULUUT, PITNESS OR AND AFFRICANAULUUT, TOTAL MARANTI, ON UNDERLES AND AFFRICANAULUUT, PITNESS OR AND AFFRICANAULUUT, PITNESS OR AND AFFRICANAULUUT, TOTAL MARANTI, ON UNDERLES AND AFFRICANAULUUT, PITNESS OR AND AFFRICANUULUUT, PITNESS OR AFFRICANUULUUT, PITNESS OR AFFRICANUULUUT, PITNESS OR AND AFFRICANUULUUT, PITNESS OR AND AFFRICANUULUUT, PITNESS OR AFFRICA

© 2013 Life Technologies Corporation. All rights reserved. The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners.

## For support visit lifetechnologies.com/support

### lifetechnologies.com

7 October 2013